


SCION PWOC (1041) Spare Parts

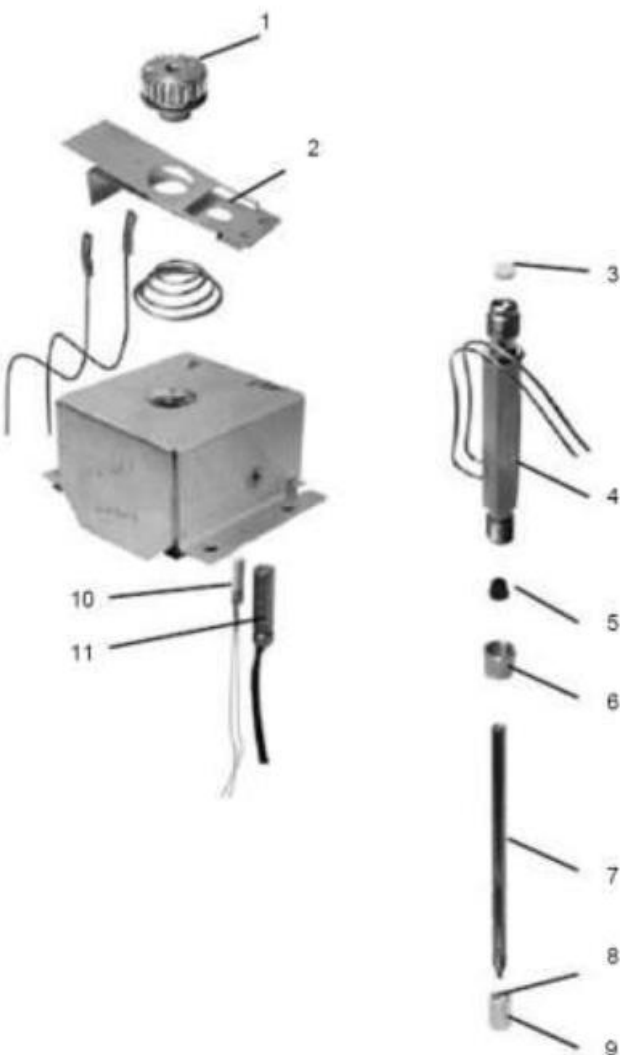

#	Description	P/N
1	Injector Nut, Autosampler	390812700
	Injector Nut, manual	392595501
	Injector Nut Wrench	390842300
2	Injector Start Switch	390820601
	3	Septum, 9.5 mm
3	BTO, lowest bleed	CR298705
	Marathon, autosampler	CR239778
	Advanced Green 3	CR246124
	Septum Pick	7200008400
3	Automatic Start Switch	390820601
4	Injector body	392548201
	5	Injector Body Ferrule
	40%G/60%V, upto400°C	CR213400
	Graphite, upto450°C	CR211400
6	Injector Body Nut, stainless steel	SWSS4021
7	0.53mmInjectorInsert	392543101
8	Column Ferrule	See ferrules
	9	Column Nut
	0.9mm Capillary, brass	394955100
10	Temperature Probe	392537401
11	Heater, 230V 125W	392539602
	Packed column Adapter kit	392558891